

Operator Settings Menu

Tactrol2[®]




A number of control system settings can be accessed via a “Hidden Menu” on the control panel, which is accessed as follows:

First turn & hold the **thermal lock key** in the override position.

Next, press the time **up** or **down** keys and release the thermal lock key.

1 should now be displayed on the temperature display and by default **0** will displayed on the time display.

Using the temperature up key scroll through from **1** to **13** in the temperature display.

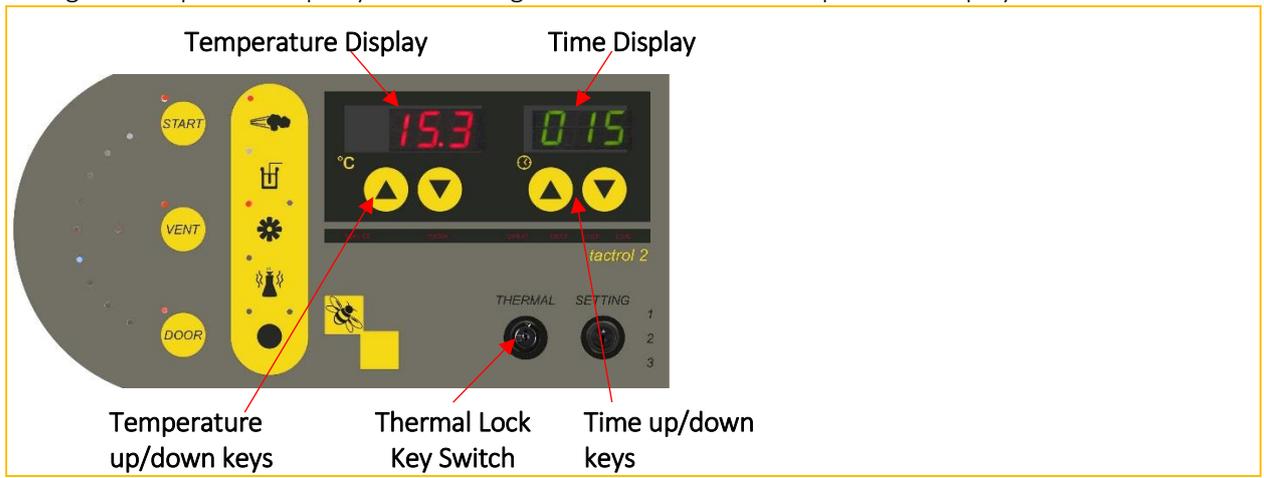

The function and setting of the individual settings are as follows:

Temp. Display	Time Display	Function	Action
Delayed Start Setting:			
01	0-24	Delayed Start Time Hour*	Enter required start time hour. (24 hour clock).
02	0-60	Delayed Start time Minute*	Enter required Start time minute.
03	0/1	Start Delay Select	1= ON, 0= OFF
Printer Settings#			
04	0-999	Print Interval (in minutes)	Enter time (minutes) between printing during process time (0= printer disabled).
Internal Clock Time Setting:			
05		Year Setting	Enter Year.
06		Month Setting	Enter Month.
07		Date Setting	Enter Day of month.
08		Hour Setting	Enter Hour. (24 Hr Clock.)
09		Minute Setting	Enter Minute.
10		Second Setting	Enter Second.

Scroll back up to parameter 1 to confirm the new or current time settings.

11	Number of Repeat Cycles [#]	Scroll up to number of repeats required
12	Service Counter Reset	Reset to zero
13	Double Door End Change [#]	Refer to separate sheet for details



After no keys are pressed for eight seconds the display returns to normal.



*For the Delay Start Timer, the time is set in real time; therefore the Main Board clock has to be correctly set for this to work properly.

After one delayed start operation, delayed start automatically switches off, and the autoclave returns to normal operation

These settings only apply where the autoclave is fitted with this functionality.