

Tactrol USB Interface Socket



The Tactrol USB interface socket is intended to allow the Autoclave user to easily download the Priorclave Tactrol Configuration (ptc) file and logged data for back-up and offering to Priorclave technical support for fault evaluation.

Using the supplied Priorclave USB Memory stick, the socket will function as follows:

Please note that the files supplied pre-installed on the memory stick must NOT be deleted, and that all files will only be used from the ROOT directory of the memory stick. It will not search subdirectories contained within the memory stick.


Other files can exist on the memory stick without causing any problems to the upload/download procedure.

Whilst the upload or download procedures are in operation it is imperative that the memory device is not removed. Removal of the device may leave the unit partially updated and may also corrupt the FAT table on the memory stick leaving it unreadable.


Creating a new USB stick for use with the available USB port:

To allow downloading Data Log Files:



1. With a new, blank Memory stick, insert it into an available PC
2. On the PC, go to the *My Computer* directory

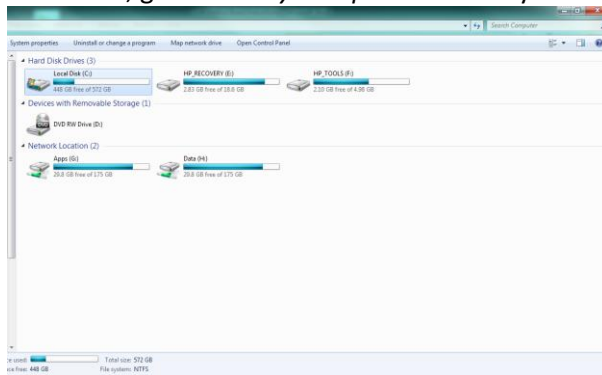


3. Double click on the appropriate memory stick option – in the example, this is Drive I

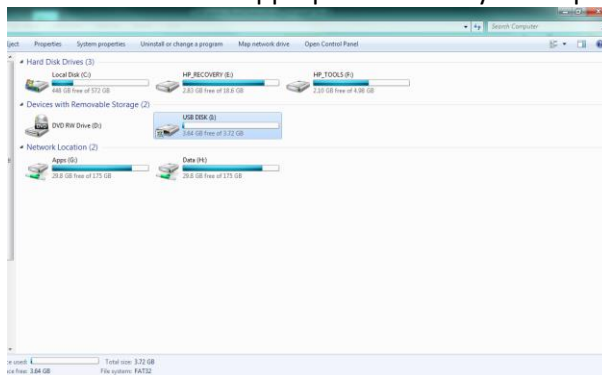


4. Open the new drive
5. Right-click on the screen and select NEW / TEXT DOCUMENT

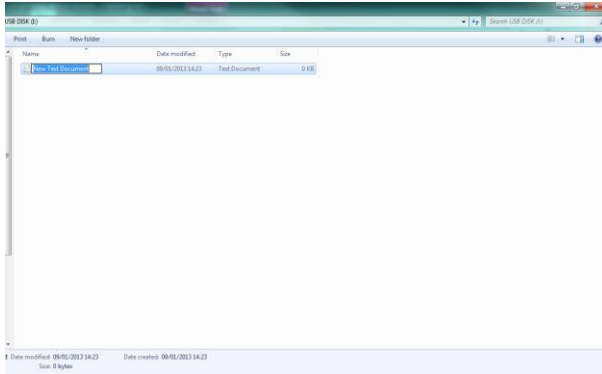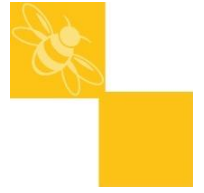


6. Name the new document 'Sxxxx' where xxxx is the Serial Number of the Autoclave

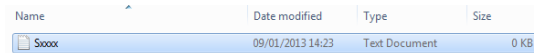

7. Open the new Text Document
8. On the first line of the document, enter the text 'DOWNLOAD=Sxxxx' where xxxx is the Serial Number of the Autoclave

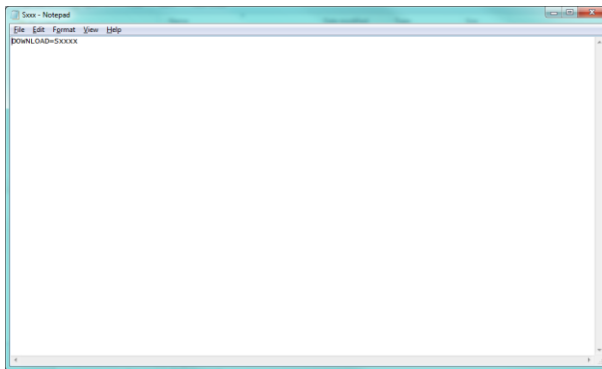

9. On the Options tab at the top of the TEXT Document window, select File / SAVE AS
10. Select the File Type option as 'All Files'
11. Save the file as name Sxxxx.DWN where xxxx is the Serial Number of the Autoclave

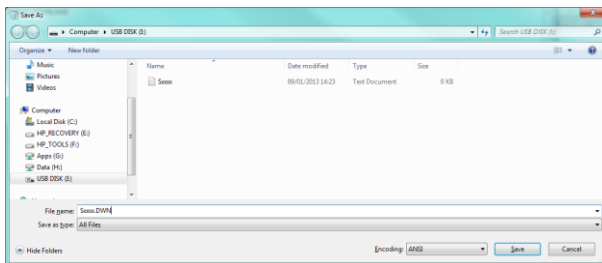

12. Press RETURN to save.

The Memory Stick is now ready to receive data log files from the Autoclave.

File transfer instructions

Retrieving a file:

To retrieve the current settings out of the Tactrol control system (ptc file), the flash drive should be attached to the USB port and power to the Autoclave must be ON.



Transfer and connection will be indicated to the user via the blue cooling LED on the main display panel. This will flash whilst the data is transferred.

Note: This operation cannot be started whilst a cycle is in progress. If the USB flash drive is plugged in whilst a program is running it will need to be removed and re-inserted when the program has completed and the unit is idle.

On successful reception of the data from the unit, the PTC file will be created and saved on to the flash drive with the file name **Sxxxx.bak**, where the serial number of the unit will be saved in the file name (marked 'xxxx').

The date and time for the file will be retrieved from the main card on-board real time clock and therefore the time and date can be checked by looking at the file time/date.

Note: Previous Sxxxx.bak files stored on the memory stick *will be overwritten*.

Updating the Autoclave Settings:

In order to update the unit settings the current setting are first retrieved as described above. This will always occur and any *Sxxxx.bak* file already in existence on the USB drive will be overwritten. The data to upload to the unit should be saved on the memory stick as a file named with the serial number of the unit to which the setting will pertain. For example, if the unit to be updated is Serial number: 1234, the file should be called S1234.ptc

Simply insert the memory stick into the socket and the new ptc file will automatically upload as indicated by the flashing blue LED. When the LED is extinguished, the uploading has been completed.

Downloading the unit's log data:

Note that, due to potential file sized, it can take several minutes to download log files.

To download the logged data the following file should be set-up on the memory stick: **Sxxxx.dwn**, where 'xxxx' represent the unit's serial number. This is originally set up as a txt file.

If the reception is successful, the following files will be created on the memory stick:

1. Sxxxx.txt (normal text file readable in any program such as notepad or imported as a tab separated file into Excel)
2. Sxxxx.enc (encoded log file that can be read by the Tactrol Log reader program or Tactrol tools log reading mechanism)

Erasing the internal log from the main card:




The internal log stored in the memory on the main card can be erased by the use of a file on the memory stick.

The following file should be present on the memory device in order to erase the log:

Sxxxx.ers

Inside this file the following text should be present: **ERASE=Sxxxx** as follows:

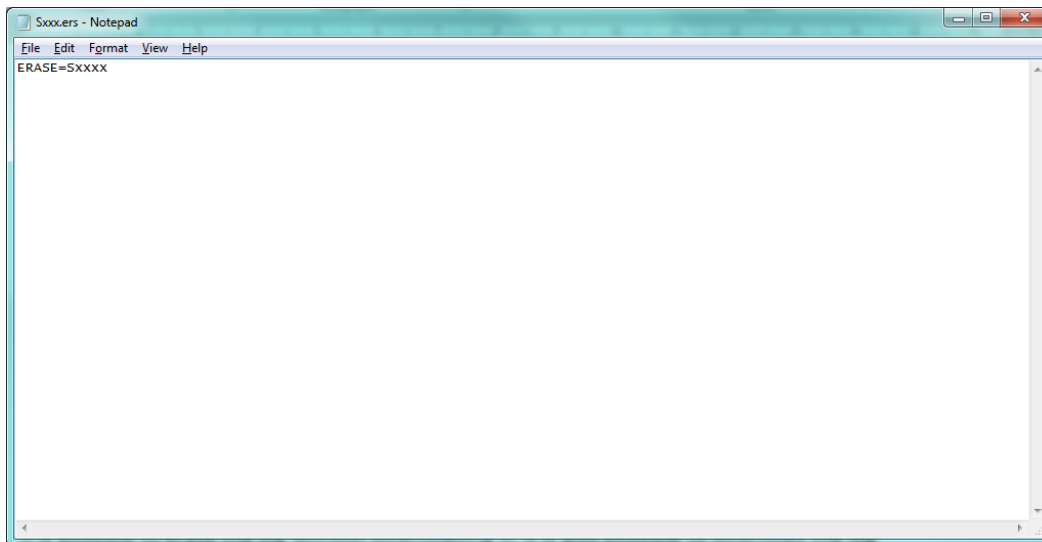

Note: no other text should be present within this file. If other text is present the erase will not be carried out.

It is possible to erase the log without downloading it; it is also possible to download the log without erasing it afterwards. However, if the log is to be downloaded then erased both files should be present. A log will always be downloaded prior to it being erased when both files are present on the memory device.

On successful erasing of the internal log, the string – LOG ERASED will be appended to the Sxxx.ers file. Therefore an erased file will look as follows:

ERASE=Sxxxx – LOG ERASED

Once this string is appended it will not erase any further logs until the – LOG ERASED text is removed from the file.