

Replacing a Paper Roll



Access to the printer for changing the paper roll is provided by a hinged front panel. The spring-loaded catch on the left-hand side of the printer front panel should be depressed to release the front panel.

The paper roll carrier is mounted on the rear of the printer front panel and will swing out from the main body for ease of roll changing.





Use the correct paper. Suitable paper is available from Priorclave Service or your local agent.

Changing Printer Ribbon

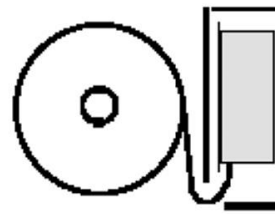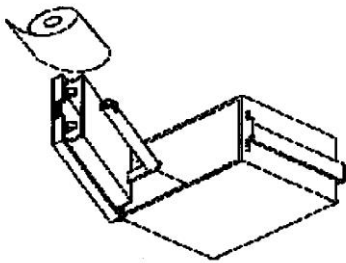

Correct Paper Path

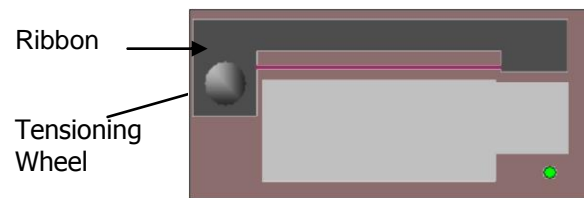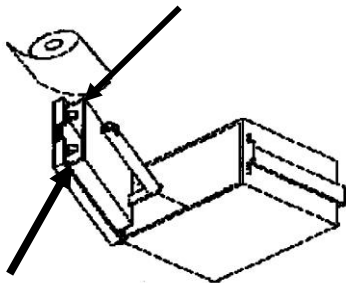

Ribbon fitting detail

The paper should pass between the ribbon and the bottom of the ribbon cassette.

The two parts of the door can now be pushed together and will click into position.