


Water Softener Manual Regeneration

If your salt unit does run out of salt you can manually regenerate the unit after adding salt and waiting 30 minutes, or you can wait for it to go through the regeneration automatically.

Firstly remove the large top cover of the Water Softener to reveal the regeneration unit:

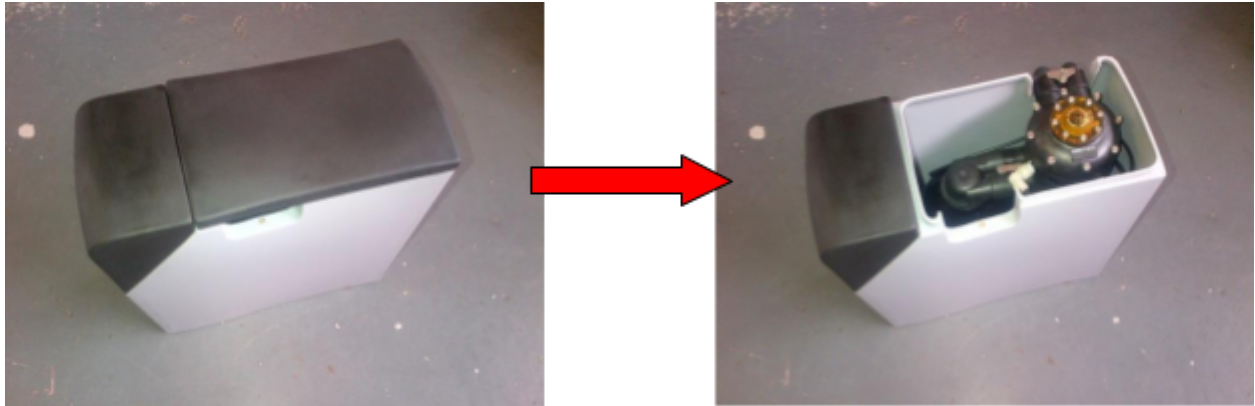

Locate adjustment screw of the regeneration unit



Using a Cross Head Screwdriver (recommended size PH2), push down firmly on the softener valve and slowly turn **CLOCKWISE** until the actuator has advanced the indicator **DOT** the 'BRINE' position. You should hear at least five clicks while turning the screw before the indicator **DOT** reaches the 'BRINE' position. At that point you will hear water flowing to the drain. This indicates that you have successfully initiated the start of the regeneration process.

Note: If you do not hear a flow of water to the drain, contact your authorized Kinetico dealer. Repeat the procedure for manual regeneration after the water flow stops (approximately 11 minutes) to be sure both resin tanks are regenerated.