



Opening Instructions

Your Priorclave Autoclave is electrically locked. In order to open it and access the full operating instructions it will first be necessary to plug it in to a suitable power supply.

After connecting the autoclave to a suitable power supply it can be opened as follows:

1. Check electricity supply is **ON**, and that the power switch is switched on.
2. When opening the autoclave for the first time the error code **F004** will be showing in the temperature display. You may ignore this at this stage as opening the autoclave and filling it with water will reset this error. Press the **door** button on the control panel there will be a bleep. Wait for about 20 seconds short time until there is another bleep and the **door** indicator illuminates. The **door** button can now be pressed again to release the lock.
3. Lift the locking handle until it reaches its safety stop. Do not lift handle until the lock has withdrawn as this may lead to damage.
4. Release the safety catch by pushing it forwards with your thumb and lift the locking handle to the top of its travel. The door can now be opened.