

Load Sense Process Timing Option

The control of the Autoclave cycle remains with the Control Probe temperature EXCEPT for the START point of the Dwell stage.

At that point, the Load Sense Process Time probe (sometimes referred to as the Wandering Probe) will start the Process Timer only when it achieves the pre-set Dwell Temperature (taking into account any offsets listed on the program settings).

While the Autoclave will control following the **RED** temperature trace, the Load Sense Process Timer probe will naturally follow a temperature trace similar to the **BLUE** line. When the Load temperature reaches that of the Dwell Set-point, marked with the 'X', the Process Timer will start.

