


## 2-Bolt RSC Door Operating Instructions

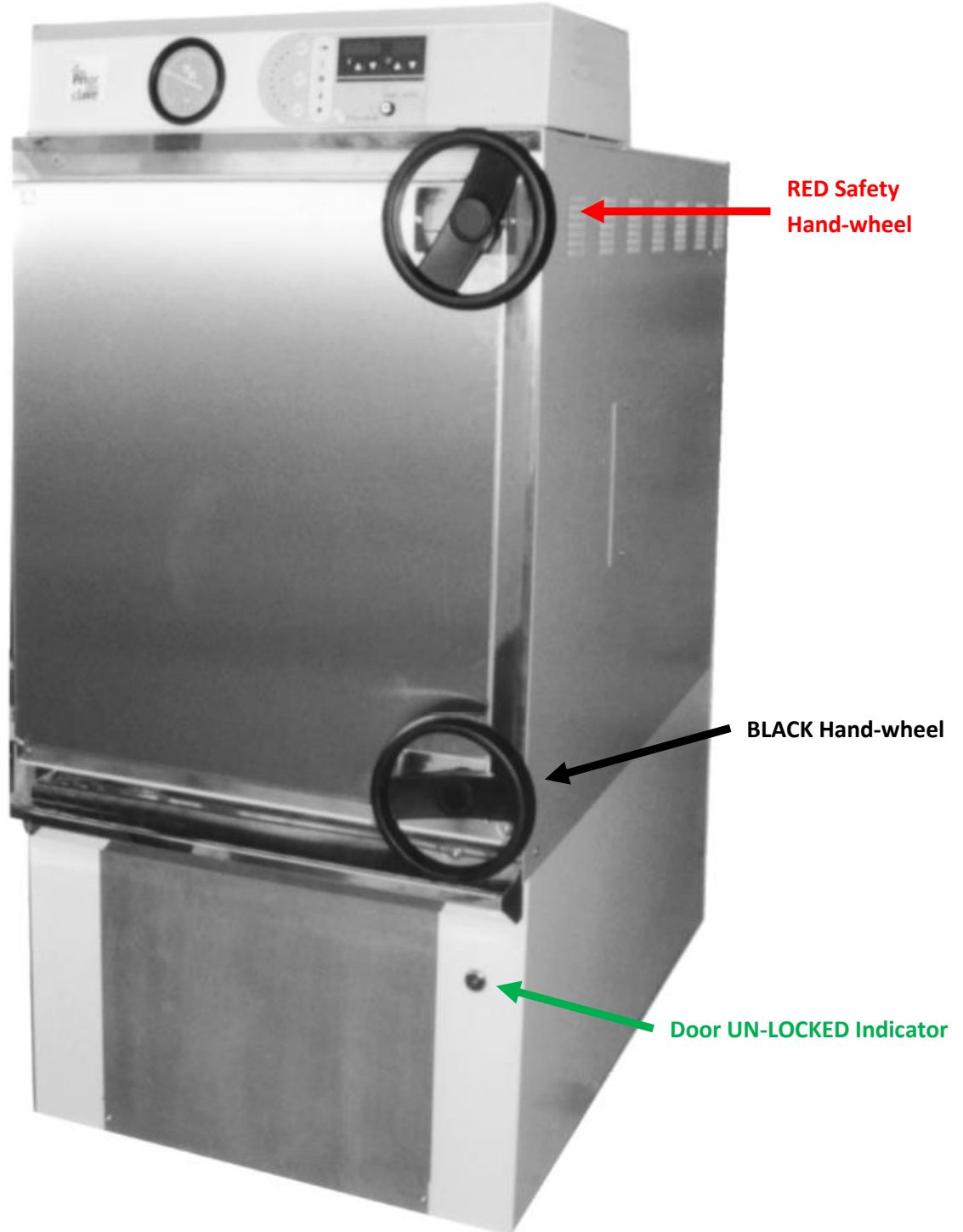


Check that the emergency stop button is not pressed in and release if necessary.

## To OPEN:

1. Check electricity and, if applicable, water supply is **ON**.

2. Press the '**DOOR**' button on the control panel there will be a bleep and the message '**hold**' will be displayed in the timer display. Wait for a short time until the timer display returns to normal, there is another bleep and the '**DOOR**' indicator illuminates.  
The '**DOOR**' button can now be pressed again to release the lock. When the lock is released a green indicator lamp below the lower hand-wheel will illuminate.  
Note that after a short time the Door Lock for the lower hand-wheel will re-engage for safety reasons.

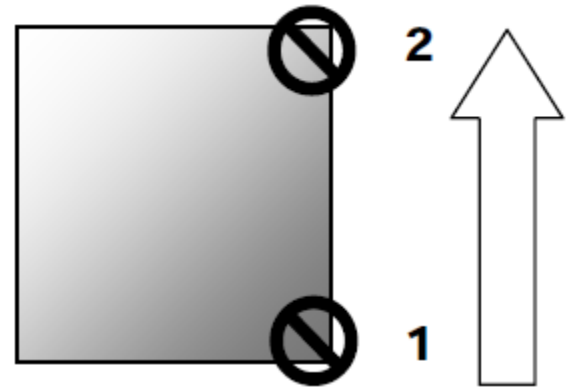

3. Open the door by undoing the lower black hand-wheel first followed by the red safety hand-wheel. With the door free, carefully pull open.


## With the Door OPEN, you can:

4. Inspect the water level - Top up with water if necessary to just below the level of the weir. If an Automatic Water-fill system is fitted, allow the autoclave to fill automatically to this level.
5. Load the autoclave directly onto the shelves or with baskets or containers.

## To CLOSE:

6. Carefully close the pressure door and start to wind in the TOP safety hand-wheel. **Tighten the wheel with hand pressure initially.**

7. Press and release the '**DOOR**' button to release the door lock. There will be a bleep and the lock will immediately withdraw. The green indicator lamp below the lower hand-wheel will illuminate whilst the lock is withdrawn. Secure by fully tightening the lower Black hand-wheel. If any difficulty is encountered, open the door and check for obstruction.  
**Tighten the wheel with hand pressure initially.**

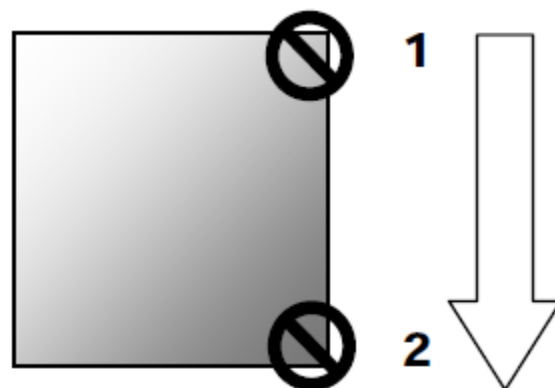

8. Once both hand-wheels are in place and hand-tight, apply one further 'nip' of the hand-wheels to fully secure.