



3-Bolt RSC Door Operating Instructions

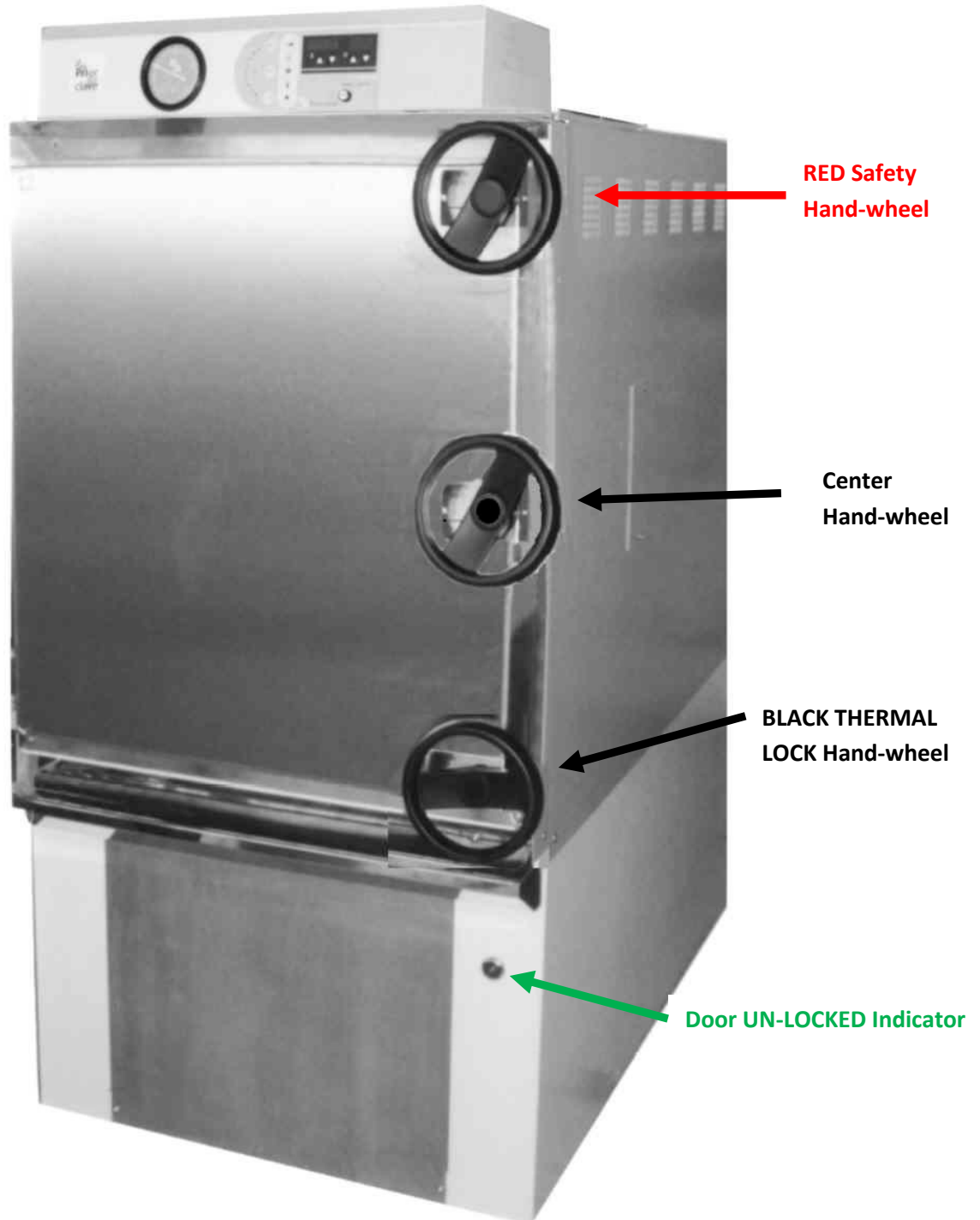

Check that the emergency stop button is not pressed in and release if necessary.

To OPEN:

1. Check electricity and, if applicable, water supply is **ON**.
2. Press the '**DOOR**' button on the control panel there will be a bleep and the message '**hold**' will be displayed in the timer display. Wait for a short time until the timer display returns to normal, there is another bleep and the '**DOOR**' indicator illuminates. The '**DOOR**' button can now be pressed again to release the lock. When the lock is released a green indicator lamp below the lower hand-wheel will illuminate.
3. Open the door by undoing the lower black hand-wheel first followed by the centre black hand-wheel and finally the red safety hand-wheel. With the door free, carefully pull open.

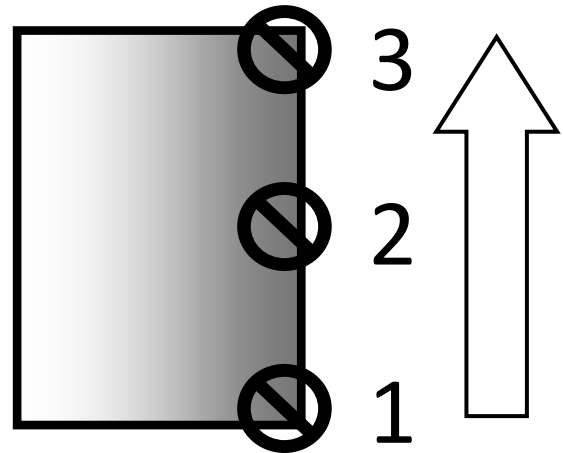


With the Door OPEN, you can:

4. Inspect the water level - Top up with water if necessary to just below the level of the weir or if an automatic. If an Automatic Water-fill system is fitted, allow the autoclave to fill automatically to this level.
5. Load the autoclave directly onto the shelves or with baskets or containers.

To CLOSE:

6. Carefully close the pressure door and start to wind in the TOP safety hand-wheel. **Tighten the wheel with hand pressure initially.**
7. Carefully close the pressure door and fully wind in the safety hand-wheel followed by the centre black hand-wheel.
8. Press and release the '**DOOR**' button to release the door lock. There will be a bleep and the lock will immediately withdraw. The green indicator lamp below the lower hand-wheel will illuminate whilst the lock is withdrawn. Secure by fully tightening the lower hand-wheel. If any difficulty is encountered, open the door and check for obstruction. **Tighten the wheel with hand pressure initially.**

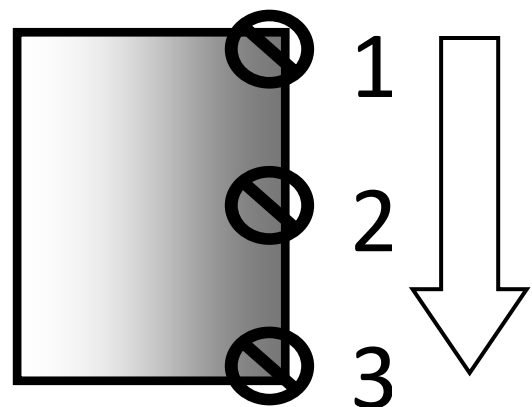

8. Once all hand-wheels are in place and hand-tight, apply one further 'nip' of the hand-wheels to fully secure.