

Tactrol Setting Key Switch Function

The Tactrol control systems can be set up to allow operate access on THREE levels using the various positions of the Setting Key Switch as follows:



Key Switch Position ONE – Key rotated fully Anti-Clockwise:

This allows only the currently selected program can be run.

Key Switch Position TWO – Key in the central position:

All programs can be selected and run. Program settings cannot be changed. Program settings cannot be changed.

Key Switch Position THREE – Key rotated fully Clock-wise:

All programs can be selected and run. Program settings can be changed freely. Program settings cannot be changed.

CANCELLING FAULT MESSAGES

Firstly, correcting the source of the original fault

Multi-program machines:

Insert the *SETTING* key into the Autoclave. Turn the key-switch to position 3, back to position 1, then back to position 3.